Germline Testing Protocol for KOMP C57BL/6N (agouti) derived ES cells (JM8A3, JM8A3.N1, JM8A1.N3)

Rev. 06/15/10

Strain Background:

- 1. The chimera's donor blastocyst C57BL/6 and the heritage is a/a (black) on the "Agouti gene" and has no Tyr^c (non-albino) on the "Color gene"
- 2. The chimera's ES cell is created from C57BL/6N with a repair at the non-agouti locus by gene targeting, and the heritage is A/a (agouti) on the 'Agouti gene', and has no Tyr^c (non-albino) on the "Color gene"
- 3. The female (C57BL/6N) breeding with the chimera is a/a (black) on the "Agouti gene" and no Tyr^c on the "Color gene"
- 4. Chimera derived from injection of the JM8 sub-line agouti (A/a) ES cells into a black host embryo (a/a) will be agouti and black. % chimerism is based on the % of agouti coat color on the mouse
- 5. These ES cells are XY and you will want to breed the male chimera. On rare occasion we encounter cells that are XØ, meaning they have lost the Y chromosome. We will notify you if this is the case for the clones or chimeras you receive. If cells are XØ you will breed the female chimera to C57BL/6N males.

Breeding chimeras with C57BL/6N females:

- 1. Set chimeras (6 wks of age), no more than one male per breeding cage, to breed with C57BL/6N females (6-8 wks of age), and no more than 3 females per cage.
- 2. If there is no sign of plug at 7 days, or pregnancy at 10 -14 days if you do not perform plug checks, remove females and give males 2-3 days rest and then replace with new females. Repeat 3 times before retiring chimera, and/or consider an artificial reproductive technique (ART) for testing chimera.
- 3. If plugs or pregnancies are observed, leave one (or both if your vivarium allows it) of the pregnant females in with male, await the birth, and allow for re-mating.
- 4. If plugs are detected but no pregnancy ensues, repeat step 1.

Germline testing:

- 1. At ten days of age, number pups (toe clip, ear tag, etc) and take tissue samples (tail snips, ear punch, etc) from **both** agouti and black pups and submit for genotyping analysis (PCR, Southern, etc).
- 2. Use germline-positive mice to establish breeding colony, and discard wild type mice. (Note: Wild type littermates should not be used for breeders as these may be host embryo derived rather than cell line derived. New C57BL/6N breeders should be requested for mating to the heterozygous positive mice for colony build, or sibling mating of heterozygous mice.)

¹ Pettitt et al.: Agouti C57BL/6N embryonic stem cells for mouse genetic resources. Nature Methods, 2009.